

2-3

Semisolid Glucose Agar.

Meat Infusion	1000 cc.	200 cc
Difco Peptone.....	10 gms.	2 gm.
NaCl.....	5 gms.	1 gm
Agar, Agar.....	2.5 gms.	5 gm.

Heat in autoclave for 1/2 hour or until the agar is dissolved.

2 gm.
Add 10 grams of dextrose. Adjust the reaction to pH 7.2 -

7.4, filter through cotton, tube and sterilize in autoclave
for 30 minutes at 15 lbs. pressure.

1